

## Exosome Isolation Kit (SEC) (0.5 mL)

JOT-EV02-03-01

### Product Overview

This product is an exosome isolation kit based on the principle of Size Exclusion Chromatography (SEC), which is a method of separating particles in solution based on particle size. The column is filled with stable and porous gel particles. When a solution containing exosomes is added to an SEC column, smaller particles are trapped in the pores, while larger molecules do not enter the pores and are eluted first. As a result, the different stages of the elution fraction will contain molecules of different sizes, starting with the larger exosome particles, followed by the smaller protein and lipid particles. This kit is more suitable for subsequent functional verification, characterization, and content detection of exosomes (nucleic acid, protein, lipid, and metabolite detection or omics analysis).

### Component and Storage Conditions

Component	Specifications	Storage Conditions
SEC Columns	0.5mL	4°C
Preservation Buffer(5x)	10mL	4°C
Balanced Buffer(10x)	10mL	4°C
Cleaning Buffer(5x)	25 mL	4°C

### Protocol

1. Open the plunger and empty the Preservation Buffer.
2. The sample loading amount is calculated based on the loading capacity of 10mg small molecule protein that can be adsorbed by 1mL of filler, and the minimum volume should not be less than 0.5mL. If the sample is less than 0.5 mL, dilute it to 0.5 mL using Balanced Buffer.
3. Balance once with Balanced Buffer (2.5 mL) that is 5 times the volume of the filler.

### QUALITY - TRACEABILITY - INTEGRITY

#### Life Science Production

The Science Centre, 1 Blaydon Road, Sandy, Bedfordshire, SG19 1RZ

Tel: +44 (0) 1234 889180 Email: sales@lifesciencegroup.co.uk Web: www.lifesciencegroup.co.uk

Life Science Production, Antibody Production Services and Diagnostic Kit Services are divisions of Life Science Group Ltd

Registered in England and Wales Company Number: 06606016




4. Load 0.5-1mL of sample, and after the sample has completely entered the filler and flowed out, add 0.5 mL of Balanced Buffer each time, collect 2-5 components with 3 drops (approximately 0.2 mL) of each component.
5. Cleaning in place: After use, clean with 10 column volumes (5 mL) of Cleaning Buffer, and then rinse 2 column volumes with distilled water (1mL), rinse 1 column volume with Preservation Buffer (0.5 mL), fill column with the Preservation Buffer, and store at 4 °C.

Usage: 20 times

## Experimental data display

Sample: Serum  
Total protein: 20mg/mL  
BCA quantification  
Sample loading: 1mL and 0.5mL  
SEC: 0.5 mL gravity column  
Collection: 1mL sample

## SDS-PAGE

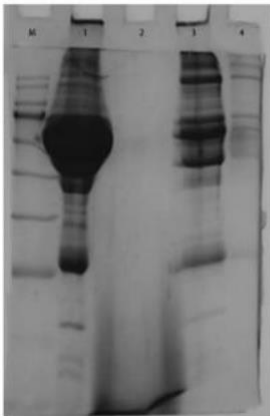

SDS: 12%  
M: 10-180KD  
Sample1: Serum 1:5 dilution  
Sample2: Cleaning in place  
Sample3: 1ml serum SEC  
Sample4: 0.5ml serum SEC

**QUALITY - TRACEABILITY - INTEGRITY**

**Life Science Production**

The Science Centre, 1 Blaydon Road, Sandy, Bedfordshire, SG19 1RZ


Tel: +44 (0) 1234 889180 Email: sales@lifesciencegroup.co.uk Web: www.lifesciencegroup.co.uk

Life Science Production, Antibody Production Services and Diagnostic Kit Services are divisions of Life Science Group Ltd

Registered in England and Wales Company Number: 06606016



## Western blot (primary antibody CD9-mAb1:2000)

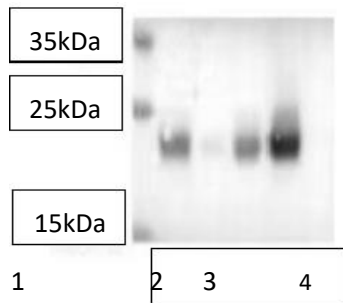

M


Sample 1: Serum control

Sample 2: Cleaning in place

Sample 3: 0.5 ml serum collection of 2-5 fractions

Sample 4: 1ml serum collection of 2-5 fractions

## ELISA test

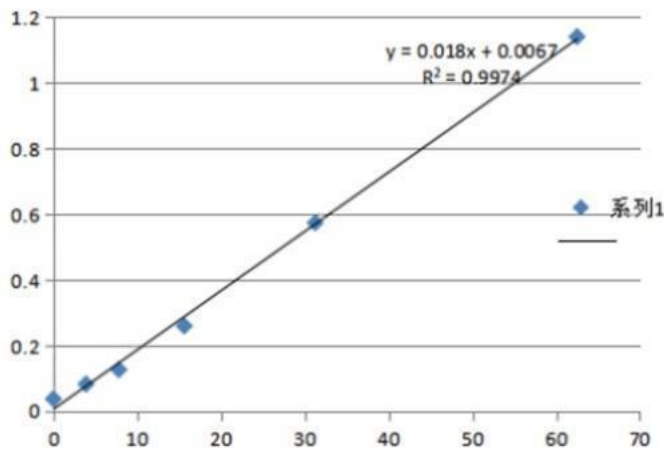

Sample	1	2	3	4	5	Balanced Buffer	Balanced Buffer
OD	0.051	0.093	0.09	0.076	0.074	0.052	0.051
Concentration	2.46	4.79	4.62	3.85	2.51	3.73	2.46

## QUALITY - TRACEABILITY - INTEGRITY

### Life Science Production

The Science Centre, 1 Blydon Road, Sandy, Bedfordshire, SG19 1RZ

Tel: +44 (0) 1234 889180 Email: sales@lifesciencegroup.co.uk Web: www.lifesciencegroup.co.uk

Life Science Production, Antibody Production Services and Diagnostic Kit Services are divisions of Life Science Group Ltd

Registered in England and Wales Company Number: 06606016



## Notes and Disclaimer

If the volume of the sample is too large after passing through the column, ultrafiltration concentration can be used to reduce the volume (Milipore ultrafiltration tube, product number: UFC801096).

Preservation Buffer (5 x), Balanced Buffer (10 x), and Cleaning Buffer (10 x) all need to be diluted to 1 x with deionized water for use.

When using, try to ensure that there is liquid inside the column to ensure that the filler can be fully wetted.

To avoid column blockage, both the buffer and sample must be treated with 0.22 µm filter membrane.

After use, the upper end of the column needs to be sealed with a sealing film to prevent the Preservation Buffer from evaporating.

Different sample volumes and chromatographic columns of different models have different ranges of component collection. Collect 1 component per well using a 96-well plate, and fluorescence intensity tests were performed to analyse the specific collection range of collected components.

This product is only for scientific research by professionals and should not be used for clinical diagnosis or treatment, nor for food or medicine.

## QUALITY - TRACEABILITY - INTEGRITY

### Life Science Production

The Science Centre, 1 Blaydon Road, Sandy, Bedfordshire, SG19 1RZ

Tel: +44 (0) 1234 889180 Email: sales@lifesciencegroup.co.uk Web: www.lifesciencegroup.co.uk

Life Science Production, Antibody Production Services and Diagnostic Kit Services are divisions of Life Science Group Ltd

Registered in England and Wales Company Number: 06606016

